



sartorius


Development of multi-spheroid co-culture 3D tumour assays using real-time live-cell analysis

K. Patel, M. Oliver, N. Holtz, T. Jackson, N. Dana, T. Dale, D. Trezise
Essen BioScience, Welwyn Garden City, AL7 3AX UK or Ann Arbor, MI, 48108, USA

Summary & Impact


- Recapitulating the tumour microenvironment (TME) is widely acknowledged as being key to the development of more predictive 3D *in vitro* efficacy models.
- Here we describe a simple & robust 3D extracellular matrix (ECM)-based workflow for culturing tumour multi-spheroids (MS) with either stromal cells or immune cells.
- IncuCyte's S3 DF-Brightfield (DF-BF) image acquisition tool enables the ability to monitor & quantify changes in spheroid size & morphology (brightfield) as well as viability (fluorescence) using real-time live-cell analysis.
- These validation methods & data demonstrate the ability to kinetically visualise & quantify immune cell-mediated toxicity within tumour multi-spheroids as well as assess the impact of stromal cells on tumour resistance to chemotherapeutic agents.
- More advanced spheroid models, incorporating the ECM & additional cell types have the potential to provide more relevant translational models for the study of the tumor micro-environment on tumour biology.

Herceptin-induced ADCC in HER2-positive multi-spheroids

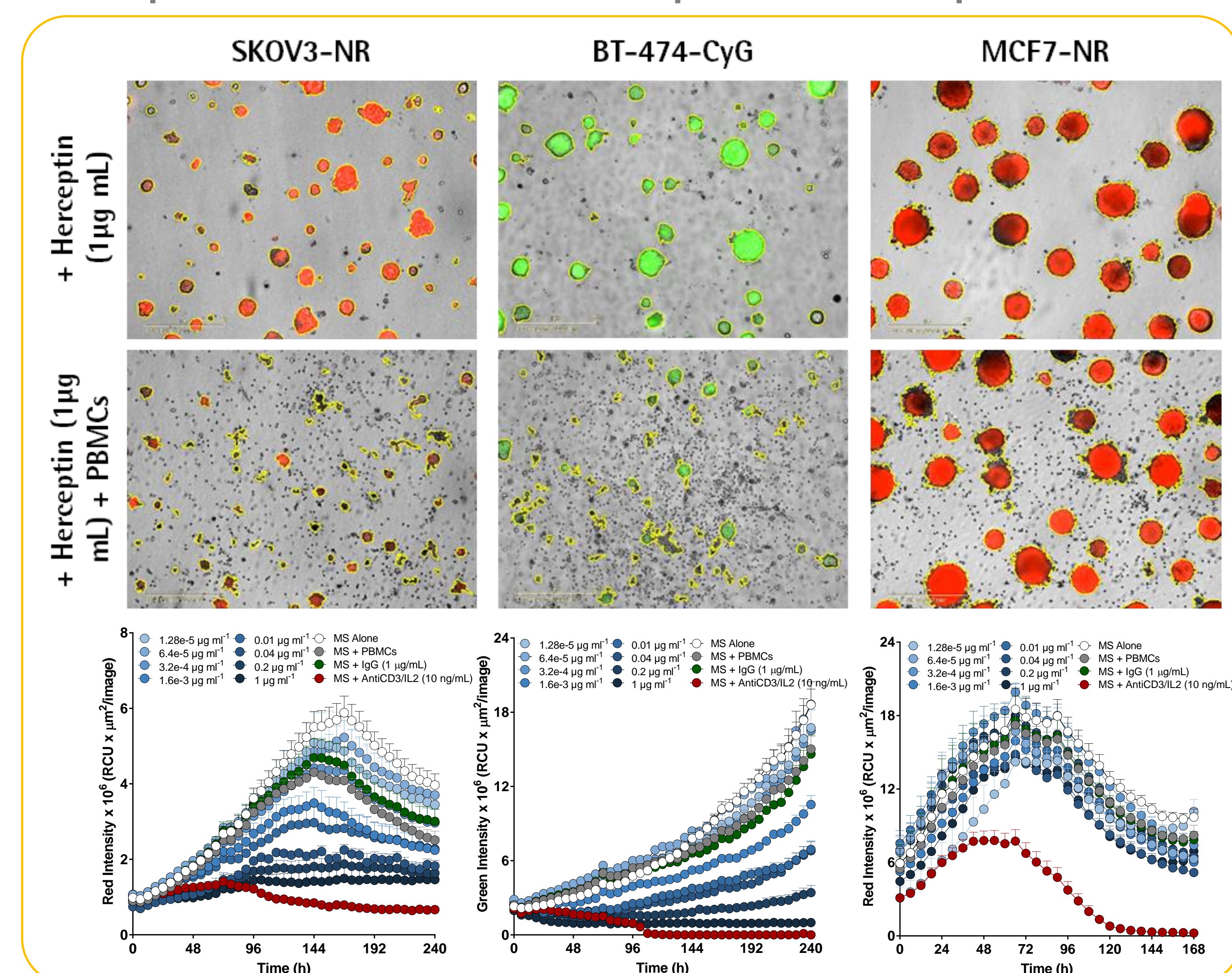

- Tumour multi-spheroids (MS) either stably expressing nuclear restricted RFP (NR) or cytoplasmic GFP (CyG) were co-cultured with PBMcs (E:T, 5:1) & treated with Herceptin (mAb targeting HER2 receptors).
- MS proliferation & immune cell-mediated cytotoxicity was quantified kinetically using the IncuCyte® S3 metrics (fluorescence intensity within BF boundary) which does not require masking of fluorescent tumour MS.
- Herceptin induced a concentration-dependent cytotoxic effect in HER2-positive (SKOV3 & BT-474) but not HER2 negative (MCF7) MS.
- BT-474 MS appear more sensitive to Herceptin cytotoxicity with the highest test concentration of Herceptin (1 $\mu\text{g ml}^{-1}$) producing ~ 80% inhibition compared with ~ 50% in SKOV3 MS.


Differential pharmacology exemplified

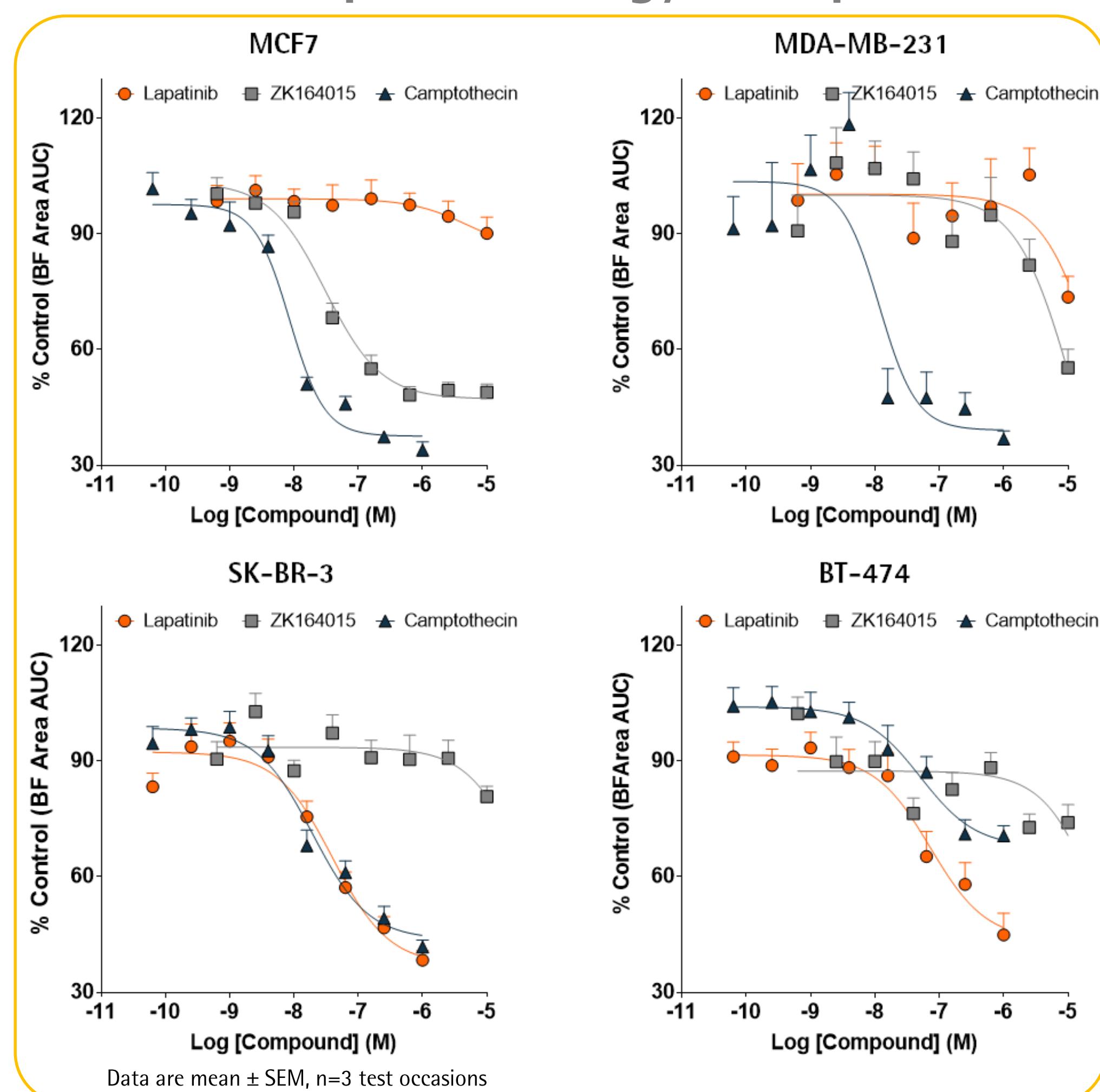


- Four tumour breast cell lines were co-cultured with normal human dermal fibroblast (NHDF's) at a 1:1 ratio (1K cells per well of each).
- MS were allowed to form for 3 days prior to treatment (7 – 10 days) with standard of care & cytotoxic agents.
- Concentration response curves represent area under curve (AUC) analysis of time-course data.
- Inhibition of MS growth caused by Lapatinib & ZK164015 aligned with the known expression profile of receptors targeted by these agents.
- The dual EGFR and HER2 tyrosine kinase inhibitor Lapatinib caused a concentration dependent inhibition of SK-BR-3 & BT-474 MS growth, while the estrogen receptor (ER) antagonist ZK164015 was a potent inhibitor of MCF7 MS growth with little or no effect on MS devoid of ER expression.
- The DNA topoisomerase inhibitor camptothecin (CMP) caused comparable inhibition of growth across all MS.

¹ Subick et al., Molecular characterization of breast cancer cell lines. Basic and Clinical Research 2010;4:35 – 41.


IncuCyte® System for continuous live-cell analysis: Methodology

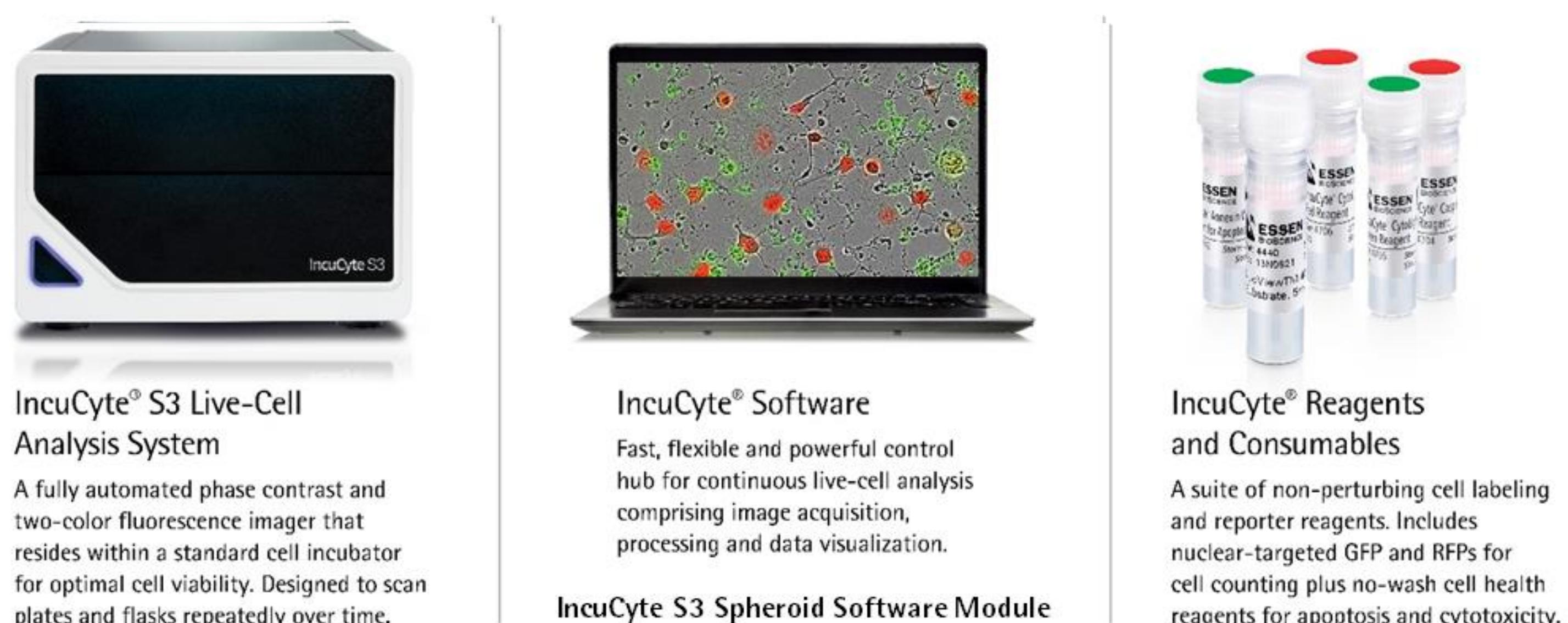


Assay workflow

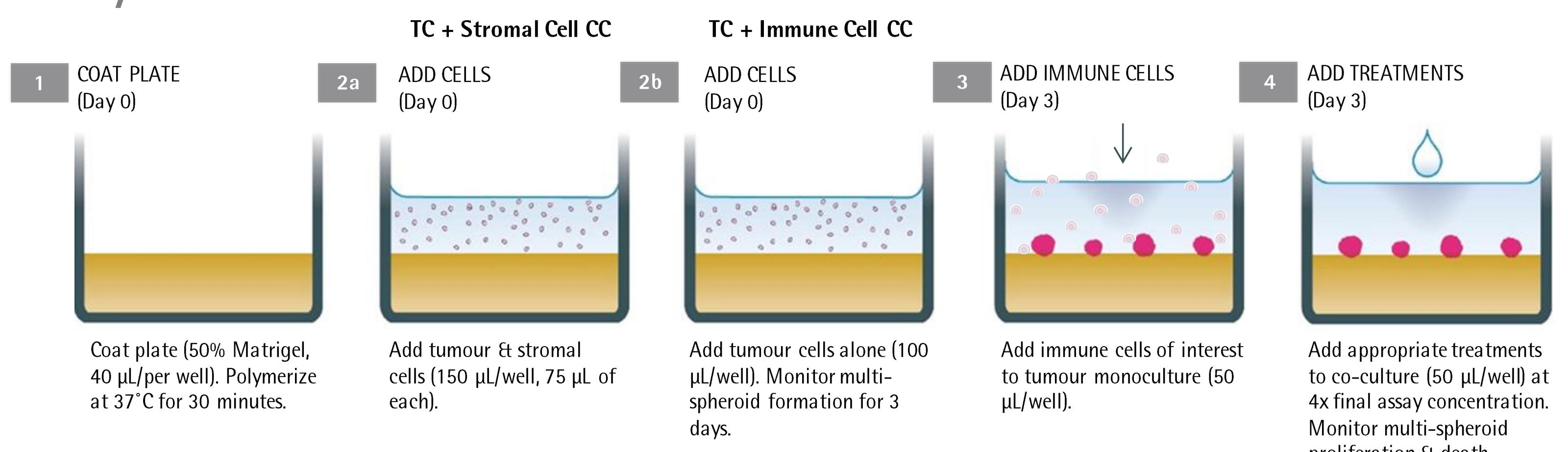

Marked effects of NHDF's on multi-spheroid morphology

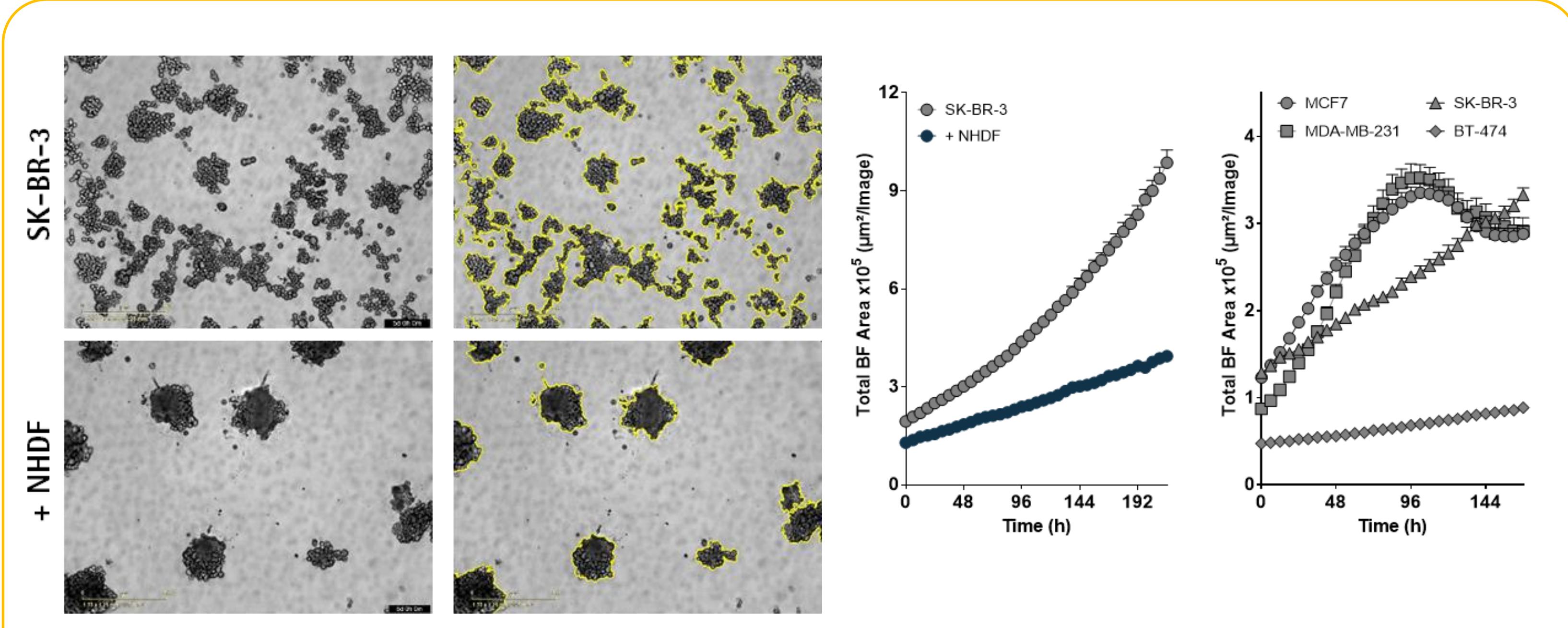


- SK-BR-3 cells were seeded alone or with NHDF's (1:1 ratio, 1K cells per well for each) & images (DF-BF) acquired every 6h. Images shown are captured 8 days post cell seeding.
- In mono-culture, SK-BR-3 cells fail to form compact multi-spheroids. Co-cultured with NHDF's results in the formation of more compact aggregates.
- IncuCyte S3 multi-spheroid software enables kinetic quantification of both morphologies, tracking MS growth via size measurements (Total BF Area).
- Note, the different growth profiles of the four tumour breast cell types in co-culture with NHDF's.

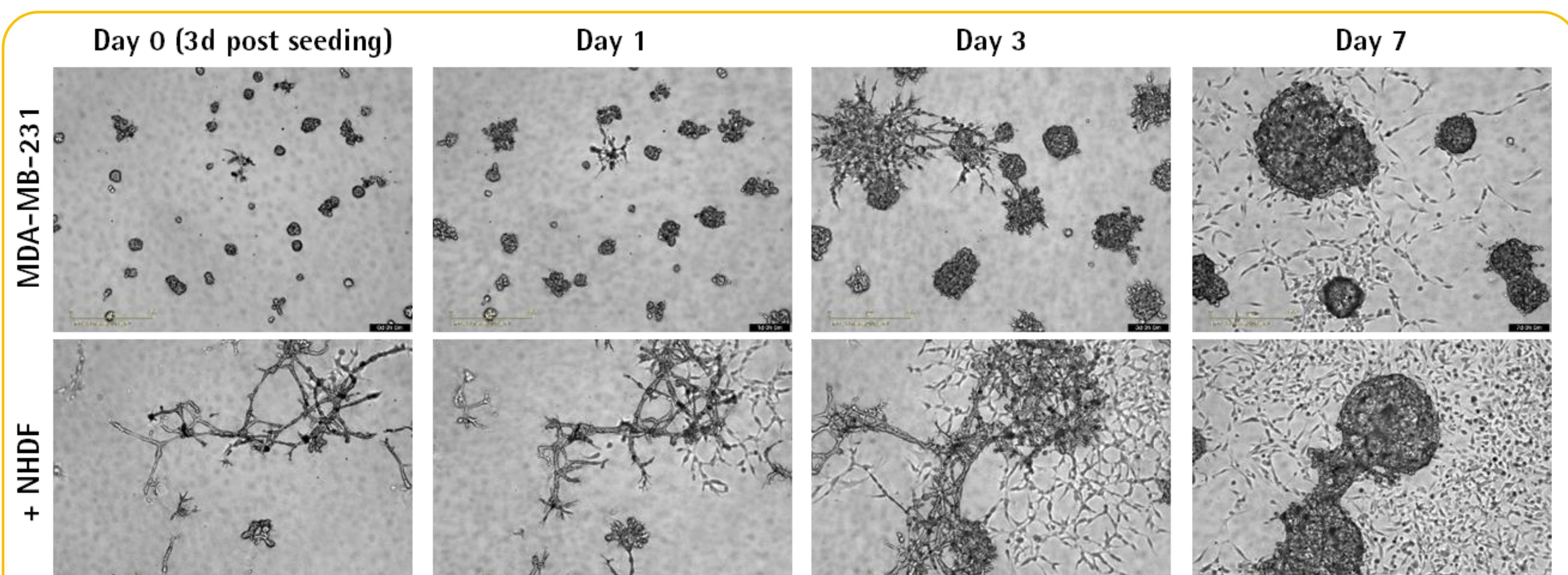

- NHDF's co-cultured with MDA-MB-231 cells (1:1, 1K per well for each) reveal a time-dependent effect on MDA-MB-231 spheroid morphology.
- MDA-MB-231 multi-spheroids transition from a stellate, branched appearance evident at early time-points, before reverting to a more clustered round appearance beyond 6 days.

Real-time quantitative assessment of treatment effects

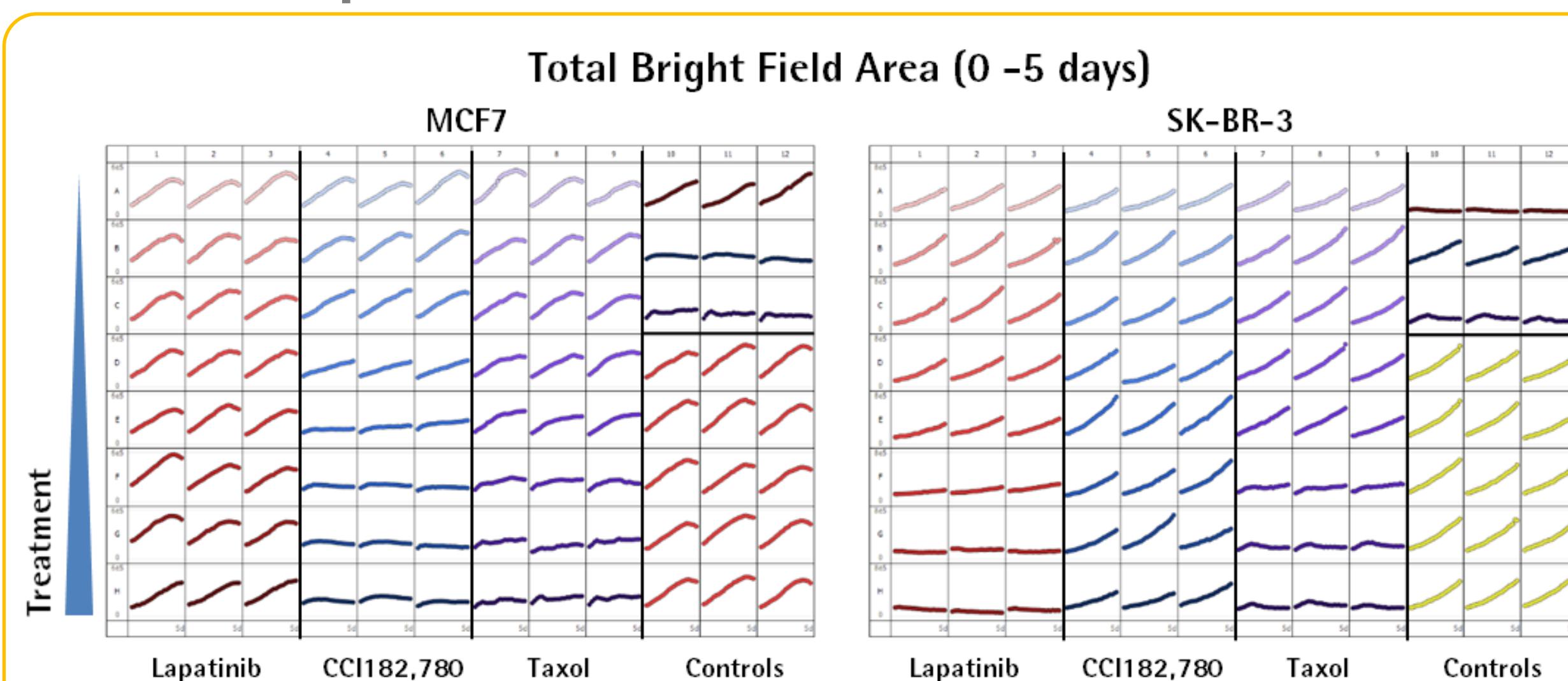

- Multi-spheroids were allowed to form for 3 days prior to treatment (5 days) with known standard of care and cytotoxic agents.
- Real-time, automated time-course plate-views enable rapid visualisation of treatment effects on MS size (Total BF Area).
- Note, the difference between cell types to treatments (time & sensitivity).