

Incucyte® Label-Free Cell Proliferation Assay

For Counting and Confluence Measurements of Adherent or Non-Adherent Cell Lines

This protocol provides an overview of the Incucyte® Label-Free Cell Proliferation Assay methodology. It is compatible with the Incucyte® Live-Cell Analysis System for kinetic, label-free analysis of cell confluence or cell counts using your choice of cells and treatments. The highly flexible assay format can be combined with our range of Incucyte® cell health and viability reagents for multiplexed measurements of cytotoxicity and apoptosis alongside proliferation in the same well.

Required Materials

- Flat bottom tissue culture plate (e.g., Corning Cat. No. 3595)

Optional Materials


- Incucyte® Cell-By-Cell Analysis Software Module (Sartorius Cat. No. 9600-0031)
 - for label-free cell counting
- Poly-L-ornithine (Sigma Cat. No. P4957)
 - for non-adherent cells
- Fibronectin (Sigma Cat. No. A7906)
 - for non-adherent cells

General Guidelines

- Following cell seeding, place plates at ambient temperature (15 minutes for adherent cell lines and 45 minutes for non-adherent cell lines) to ensure homogenous cell settling.
- Remove bubbles from all wells by gently squeezing a wash bottle containing 70-100% ethanol with the inner straw removed, to blow vapor over the surface of each well.
- After placing the plate in the Incucyte® Live-Cell Analysis System, allow the plate to warm to 37° C for 30 minutes prior to scanning.


Quick Guide

1. Coat wells

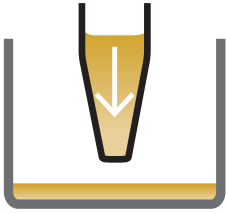


Coat wells of plate (50 μ L/well) with appropriate matrix. Optional for adherent cell lines.

2. Plate cells

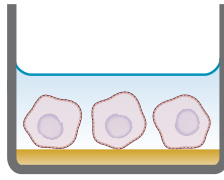

Seed cells (100 μ L/well, 1,000–10,000 for adherent and 5,000–50,000 for non-adherent) into a 96-well plate.

3. Add treatments

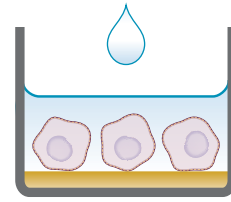

Add desired treatments (100 μ L/well, 1X for adherent cell lines, 2X for non-adherent cell lines).

Adherent Cell Line Protocol

Day 0

1. Coat Wells (optional)

- 1.1 Depending on cell line used, coat a 96-well flat bottom plate with relevant coating matrix according to manufacturer's recommendation.
- 1.2 Prior to cell seeding, prepare cell treatments at 2x final assay concentration in enough cell culture medium to achieve a volume of 100 μ L per well.

2. Plate Cells

- 2.1 Seed your choice of cells (100 μ L per well) at an appropriate density into a 96-well plate, such that by day 1 the cell confluence is approximately 10–20%. The seeding density will need to be optimized for the cell line used; however, we have found that 1,000 to 2,500 cells per well (10,000 to 25,000 cells/mL seeding stock) are reasonable starting points.
 - a. Monitor cell growth using the Incucyte® Live-Cell Analysis System to capture phase contrast images every two hours and analyze using the integrated confluence algorithm.

- 3.2 Remove the cell plate from the incubator and aspirate medium from wells.
- 3.3 Add treatments and controls to appropriate wells of the 96-well plate.
- 3.4 Place the cell plate into the Incucyte® Live-Cell Analysis System and allow the plate to warm to 37° C for 30 minutes prior to scanning.
 - a. Objective: 4X, 10X or 20X
 - b. Channel selection: Phase Contrast (+ Fluorescence if fluorescent label or cell health reagents are used)
 - c. Scan type: Standard
 - d. Scan interval: Typically, every 1 to 2 hours, until your experiment is complete.

Note: Label free cell counting can be enabled on Incucyte® Live-Cell Analysis System with use of the Incucyte® Cell-by-Cell Analysis Software Module.

- a. Scan type: Standard | Adherent Cell-by-Cell
- b. Objective: 10X

For further details of this analysis module and its application see: www.essenbioscience.com/cell-by-cell

Day 1

3. Add Treatments

- 3.1 Prepare 1X concentrations of desired cell treatments in cell culture medium. The volumes may be varied; however, we recommend preparing enough volume of each desired treatment | dilution in order to achieve 100 μ L per well.

Non-Adherent Cell Line Protocol

Day 1

1. Seed Cells and Add Prepared Treatments

- 1.1 Coat a 96-well flat bottom plate with relevant coating matrix. We recommend coating with 50 µL of either 0.01% poly-L-ornithine solution or 5 µg/mL fibronectin diluted in 0.1% BSA. Coat plates for 1 hour at ambient temperature, remove solution from wells, then allow plates to dry for 30–60 minutes prior to cell addition.
- 1.2 Prior to cell seeding, prepare cell treatments at 2X final assay concentration in enough cell culture medium to achieve a volume of 100 µL per well.

2. Plate Cells

- 2.1 Seed your choice of cells (100 µL per well) at an appropriate density into a 96-well plate. The seeding density will need to be optimized for the cell line used; however, we have found that 5,000–50,000 cells per well (50,000–500,000 cells/mL seeding stock) are reasonable starting points.

Note: If studying immune cell clustering and proliferation, prepare cell activation treatments at 5X final concentration, and immediately add 50 µL per well containing cells. It is advised that some control wells containing only vehicle are included in the plate.

3. Add treatments

- 3.1 Immediately after cell seeding, add treatments and controls to appropriate wells of the 96-well plate containing cells. Triturate wells to appropriately mix the treatment to ensure cell exposure at 1X.
- 3.2 Place the cell plate into the Incucyte® Live-Cell Analysis System and allow the plate to warm to 37° C for 30 minutes prior to scanning.
 - a. Objective: 4X (recommended 1 image per well or whole well) or 10X
 - b. Channel selection: Phase Contrast (+ “Green” or “Red” if fluorescent label or cell health reagents are used)
 - c. Scan type: Standard
 - d. Scan interval: Typically, every 1 to 2 hours, until your experiment is complete.

Note: Label-free cell counting can be enabled on Incucyte® Live-Cell Analysis System with use of the Incucyte® Cell-by-Cell Analysis Software Module.

- a. Scan type: Standard | Non-Adherent Cell-by-Cell
- b. Objective: 20X

For further details of this analysis module and its application see: www.essenbioscience.com/cell-by-cell

Related Products and Applications

A comprehensive range of fluorescent nuclear labeling and cell health reagents are available for use with the Incucyte® Live-Cell Analysis System to enable multiplexed measurements of cytotoxicity and proliferation alongside apoptosis.

Compatible with the Incucyte® Live-Cell Analysis System

Proliferation & Cell Cycle				
Product	Color	Quantity	Compatibility	Cat. No.
Incucyte® Nuclight Green Lentivirus (puro)		One vial: 0.2 mL	SX1, S3, SX5	4624
		One vial: 0.6 mL		4475
Incucyte® Nuclight Green Lentivirus (bleo)		One vial: 0.2 mL	SX1, S3, SX5	4626
		One vial: 0.6 mL		4477
Incucyte® Nuclight Red Lentivirus (puro)		One vial: 0.2 mL	SX1, S3, SX5 (Green Red Optical Module)	4625
		One vial: 0.6 mL		4476
Incucyte® Nuclight Red Lentivirus (bleo)		One vial: 0.2 mL	SX1, S3, SX5 (Green Red Optical Module)	4627
		One vial: 0.6 mL		4478
Incucyte® Nuclight Orange Lentivirus (puro)		One vial: 0.2 mL	S3 for Neuroscience, SX5	4771
Incucyte® Nuclight NIR Lentivirus (puro)		One vial: 0.2 mL	S3 for Neuroscience, SX5	4805
Incucyte® Nuclight Rapid Red Dye		One vial: 50 µL	SX1, S3, SX5 (Green Red Optical Module)	4717
Incucyte® Nuclight Rapid NIR Dye		One vial: 50 µL	S3 for Neuroscience, SX5	4804
Incucyte® Cell Cycle Green Red Lentivirus (puro)		One vial: 0.6 mL	SX1, S3, SX5 (Green Red Optical Module)	4779
Incucyte® Cell Cycle Green Orange Lentivirus (puro)		One vial: 0.6 mL	SX5	4809

Apoptosis & Cytotoxicity				
Product	Color	Quantity	Compatibility	Cat. No.
Incucyte® Annexin V Green Dye		One vial: 100 tests	SX1, S3, SX5	4642
Incucyte® Annexin V Red Dye		One vial: 100 tests	SX1, S3, SX5 (Green Red Optical Module)	4641
Incucyte® Annexin V NIR Dye		One vial: 100 tests	S3 for Neuroscience, SX5	4768
Incucyte® Annexin V Orange Dye		One vial: 100 tests	S3 for Neuroscience, SX5	4759
Incucyte® Caspase-3/7 Green Dye		One vial: 20 µL	SX1, S3, SX5	4440
Incucyte® Caspase-3/7 Red Dye		One vial: 20 µL	SX1, S3, SX5 (Green Red Optical Module)	4704
Incucyte® Cytotox Green Dye		Five vials: 5 µL each	SX1, S3, SX5	4632
Incucyte® Cytotox Red Dye		Five vials: 5 µL each	SX1, S3, SX5 (Green Red Optical Module)	4633

*Pre-labeled Nuclight cell lines are also available for purchase. Please visit shop.essenbioscience.com for more information

A complete suite of cell health applications is available to fit your experimental needs.
Find more information at www.sartorius.com/incucyte

For additional product or technical information, please email us at AskAScientist@sartorius.com

For Research Use Only. Not For Therapeutic or Diagnostic Use.

Sales and Service Contacts

For further contacts, visit
www.sartorius.com

Essen BioScience, A Sartorius Company

www.sartorius.com/incucyte

E-Mail: AskAScientist@sartorius.com

North America

Essen BioScience Inc.
300 West Morgan Road
Ann Arbor, Michigan, 48108
USA
Telephone +1 734 769 1600
E-Mail: orders.US07@sartorius.com

Europe

Essen BioScience Ltd.
Units 2 & 3 The Quadrant
Newark Close
Royston Hertfordshire
SG8 5HL
United Kingdom
Telephone +44 1763 227400
E-Mail:
euorders.UK03@sartorius.com

APAC

Essen BioScience K.K.
4th Floor Daiwa Shinagawa North
Bldg.
1-8-11 Kita-Shinagawa
Shinagawa-ku, Tokyo
140-0001
Japan
Telephone: +81 3 6478 5202
E-Mail: orders.US07@sartorius.com

Specifications subject to change without notice.

© 2020. All rights reserved. Incucyte, Essen BioScience, and all names of Essen BioScience products are registered trademarks and the property of Essen BioScience unless otherwise specified. Essen BioScience is a Sartorius Company. Publication No.: 8000-0395-H00
Status: 08 | 2020